



www.insize.com




**5309-DM51
DIGITAL MICROSCOPE
OPERATIONAL MANUAL**



1. Packing List




1

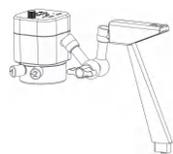


2



3

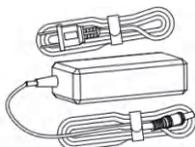

4

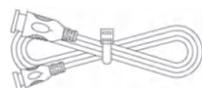

5



6



7



8



9

1. Tripod
2. Camera
3. Display screen
4. AC/DC adapter
5. HDMI cable

6. Screw
7. Screwdriver
8. Inner hexagon spanner
9. Remote controller

2. Installation Guide

2.1 Display assembly

- ◆ Unscrew one screw on the support column, as shown in Figure 1.
- ◆ Take out the tripod from the box and lock the support column on the tripod with screws and washers, as shown in Figure 2.
- ◆ Take out the display screen from the box and fix the display cover on the connecting rod with four M3 screws, as shown in Figure 3.

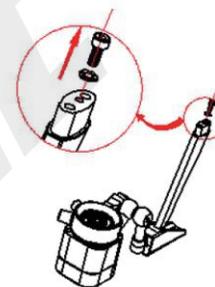


Figure 1

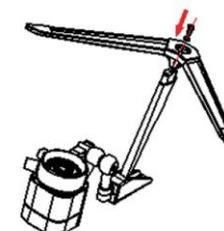


Figure 2


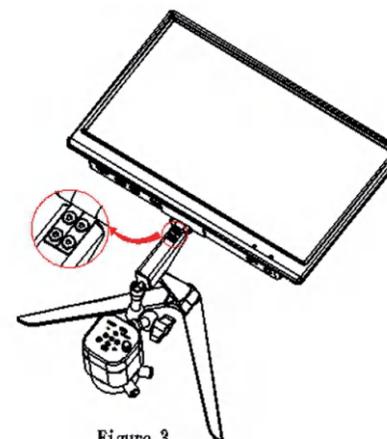


Figure 3

2.2 Cable connections

- ◆ Connect mini port of the HDMI cable with the HDMI Mini port of the camera, and fasten it with the screws provided, as shown in Figure 4; the other end HDMI-A is connected to the display monitor, as shown in Figure 4.
- ◆ Power the display screen with provided 110-240V AC plug, as shown in Figure 5.

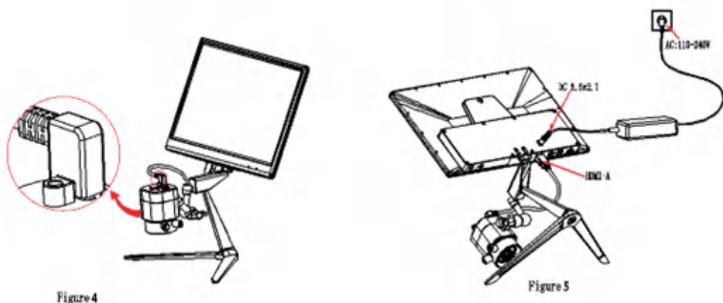

2.3 Start to use

- ◆ Turn on the monitor and camera
- ◆ Adjust the camera position
- ◆ Adjust the focal length

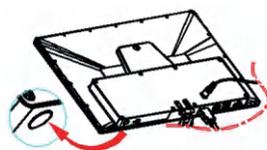


Figure 6: Turn on the monitor

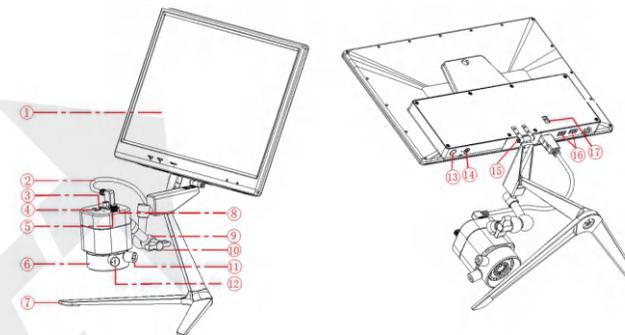


Figure 7: Turn on the camera



Figure 8: Adjust the focal length on the top of camera to produce clear image.

3. Function Introduction of Each Part



- | | |
|-------------------|---|
| 1. Display screen | 15.6" display screen, for image display and USB power supply interface. |
| 2. HDMI cable | Power supply and signal transmission between camera and display screen. |
| 3. Push-button | 5 buttons can match to each other to achieve different functions.

Left button: Menu/return/frozen (hot key) function.

Bottom button: Zoom out/downward/video fast forward function.

Top button: Zoom in/upward/speed retreat function.


Middle button: Record/camera/play back function.

Right button: Mode/sure/cross hair (hot key) function. |


- 4. digital camera Switch Put the switch to right: turn on the digital camera; put the switch to left: turn off the digital camera.
- 5. Focus knob Rotate the knob to implement the focus function; Rotate clockwise:back focal length decreases and object distance increases; Rotate anticlockwise:back focal length increases and object distance decreases.
- 6. light lamp The coaxial light and the diffuse reflection coexist, and the two effect lights are respectively controllable;thecoaxial light lamp adopts a circular distribution of 13 light beads at a small angle;the diffuse reflection light is distributed by 18 high-brightness lightbeads,and the brightness is uniform.
- 7. Tripod Aluminium alloy tripod.
- 8. TF card slot The location of TF card, photos and videos will be stored in TF card.
- 9. Stand can rotate by 360 degree.
- 10. Mobile knob Loosen the knob anticlockwise and the stand will be adjusted in any degree;tighten the knob clockwise and the current position of the electronic eyepiece can be fixed.
- 11. Adjustment knob of the coaxial light Rotate clockwise: turn on the coaxial light/strengthen coaxial lightness;rotate counterclockwise: reduce the coaxial lightness/turn off the coaxial light.
- 12. Adjustment knob of the diffuse reflectance light Rotate clockwise: turn on the diffuse reflectance light/strengthen the diffuse reflectance lightness; rotate counterclockwise:weaken the diffuse reflectance lightness/turn off the diffuse reflectance light.
- 13. Monitor switch Press and hold the on key for more than 4seconds to turn on or off.(the digital camera stops working after the monitor is turned off)
- 14. Remote controller interface Connect to the remote controller to adjust parameters of the monitor.

- 15. Rotating shaft It is a 5V 2A power output interface, provide power to digital camera through USB cable.
- 16. USB power supply interface Power supply to the monitor.
- 17. Power interface Satisfies the display up and down flip.

4. Accessory Introduction

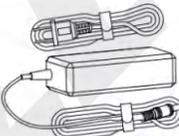- 

AC/DC adapter Connect and supply power to the monitor.
- 

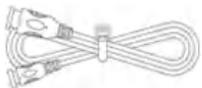Remote controller Connect with electronic eyepiece or monitor for adjusting parameters.
- 

HDMI cable Connect the electron microscope to the display can realize signal.

5. Instructions

5.1 Introduction of Various Functions

- ◆ The machine contains altogether 5 press buttons.
- ◆ Short press refers to press time 0.2s~1s. Long press refers to press time more than 2s or above. (Without special reference, press below all means short press).
- ◆ For 10 seconds without operation, the system will automatically hide the icon on the screen desk for users to watch the video more conveniently. The icon on the desk can be awakened by pressing the left button and the other 4 buttons will execute the command of the current press button while been awakened. (Users can press arbitrary button to awake the desk icon).
- ◆ The maximum magnification ratio of the zoom in function button: x4. Zoom out function button can only be available when being zoomed in first.

5.2 Photography

- ◆ The icon  on the top right corner of the screen displays for photography mode. Please ensure the TF card insertion, press “middle button” for picture shooting. Pictures will be automatically saved into the Photo file under TF card file Microscope. File name format: "Photo_00001.jpg", "Photo_00002.jpg", ...

5.3 Record

- ◆ The icon  on the top right corner of the screen displays for recording mode to record sound and video files. Please ensure the TF card insertion, press “middle button” for recording and press “middle button” again to exit recording. Records will be automatically saved into the Video file under TF card file Microscope. File name format "Video_0001.mov", "Video_0002.mov" ...
Note: The maximum capacity of recording file is 4G. When 4G capacity is reached, a new recording file will be started.

5.4 Playback

- ◆ The top right corner icon  displays for Playback mode. Flip over for observation by pressing “Top Button” or “Bottom button”. Pressing button “Left button” can enter the file attributes interface, which includes three operation functions such as delete, protect and slide show.

5.5 Frozen

- ◆ Under the mode of camera or recording, long press “left button” and the right side of the desk will display the icon  which is the frozen screen. You can long press “the left button” or short press any button to defrost the screen and the icon  disappears. Each button's function is performed after unfreeze.

5.6 Menu Setup

- ◆ Pressing “Left Button” to enter the menu bar interface. Through button “Top Button” and “Bottom Button” to move the cursor up and down for selecting the required function. Press “Right Button” to enter into setup. After setup, press “Left Button” to exit menu bar.

Note: common functions are set up on the menu bar, namely exposure compensation, white balancing, 3 sequence, image quality, HDR, resolution, sharpness, shooting ways, color, quick preview, button press sound, auto power off, language, frequency, image rotation, frozen and date/time.

5.7 Exposure setup

- ◆ Under the mode of camera or recording, press the “left button” to enter the menu bar. Press the “top button” or the “bottom button” to select the menu bar “exposure EV” and then press the “right button” to enter the exposure setup interface.

Then the screen shows the icon . (Note: font flashes: stands for the cursor is located in current position. Red font: stands for the status of been selected.)

Press the “top button” or “bottom button” to move the cursor. Press the “right button” to select the current status and the font color will turn red. Press the “left button” to exit the exposure setup interface.

The cursor points to A, the real-time automatic exposure function will be started by pressing the “right button” and the color of the letter A will turn red.

The cursor points to S, the single real-time automatic exposure function will be started by pressing the “right button” and the color of the letter S will turn red.

- ◆ The cursor points to M, the manual exposure can be started by pressing the “right button” and the color of the letter M will turn red and a flashing red number 0 will also appear on the right side. Press the “top button” or “bottom button” to change the current number value (the valid adjustment range is -10~+10 and the default value is 0). Press the “right button” to save the number value and the number value will be covered simultaneously.

5. 8 White balance adjustment

- ◆ Under the mode of camera or recording, press the “left button” to enter the menu bar. Press the “top button” or the “bottom button” to select the menu bar “white balance WB” and then press the “right button” to enter the interface of the white balance setup. Now the screen will display the icon . (Note: font flashes: stands for the cursor is located in current position. Red font: stands for the status of been selected.) Press the “top button” or the “bottom button” to move the cursor. Press the “right button” to select the current status and the font color will turn red. Press the “left button” to exit the White balance setup interface.

The cursor points to A, the real-time automatic white balance function will be started by pressing the “right button” and the color of the letter A will turn red.

The cursor points to S, the single automatic white balance function will be started by pressing the “right button” and the color of the letter S will turn red.

The cursor points to M, the manual white balance can be started by pressing the “right button” and the color of the letter M will turn red and a group of value  will also appear on the right side. The flashing red number is the value currently been selected. Press the “top button” or “bottom button” to change the current number value (the valid adjustment range is 1-255). Press the “right button” to switch the value R, G and B consequently.

Whichever value letter (R / G / B) is selected, as long as the “left button” is pressed, RGB current number value will be saved and altogether RGB and the number value will be covered.

Note: From automatic white balance to manual white balance: the value of manual white balance is the white balance under the current background. From single automatic white balance to manual white balance: the value of manual white balance is the previous single automatic white balance.

5.9 Crosshair mark

Under the mode of camera or recording, long press the “right button” to recall the crosshair mark and then long press the “right button” to cancel the crosshair mark. The default cross line invoked by shortcut key is the first kind. If either of the other two shortcut keys of the cross line is needed, other cross lines will be firstly invoked via menu and current cross line is controlled by shortcut key.

There are three kinds of crosshair marks (“crosshair mark” menu can be found in the menu bar):

1. Cross (default)
2. Grid
3. Concentric circles

Notes and Safety Requirements

Cautions and Notes

- ◆ To avoid danger or damage incurred to the lens, do not touch the lens or sensor directly with your fingers.
- ◆ To avoid failure or electric shock hazard and so on, do not disassemble or modify the internal structure of the device.
- ◆ Do not plug in or unplug the Power connection port when hands are wet.
- ◆ Do not use alcohol and other organic solvents to clean.
- ◆ If the lens or sensor is dirty or damp, you should better use dry and non-linen fabric or professional lens tissue to wipe them. To avoid scratches on the surface, do not touch the lens with your fingers. Wipe the lens or sensor lightly.
- ◆ The products are not specifically designed for an outdoor use. Do not expose it to outdoor environment without any protection. Excessive temperature and humidity will damage the lens. Please avoid using the product under the following environment: high temperature or high humidity environment, places with direct sunlight, dirt or vibration and places near heat source.
- ◆ Please use and store in the following environment:
 Operating temperature : -10°C~ 40°C
 Storage temperature: -10°C~ 60°C
 Operating Humidity: 30~60%RH
 Storage Humidity: 10~80%RH
- ◆ If any foreign matter, water or liquid enter into the device by accident, disconnect the Power line immediately. Please send it to the maintenance center and do not use the hair dryer to dry it by yourself.
- ◆ To prevent microscope from being tripped over or dropped, please put away the device's connect cable in use or standby.
- ◆ To avoid electric shock by accident, please power off microscope before you move your display or equipment.
- ◆ The cleanliness of the device lens will directly affect clarity degree of contents from the computer screen during preview. Problems like various circles or spots on the screen may mostly be incurred by dirt on the lens. When cleaning, please use professional lens tissue or other professional detergent to clear the dirt on the lens.

Disclaimer

- ◆ In order to protect the legitimate rights and interests of users, please carefully read the instructions, disclaimers and safety instructions provided with this product before using it. The company reserves the right to update the above documents. Please operate the product according to the instructions and safety instructions.
- ◆ Once you begin to use the product, you shall be deemed to have read, understood, recognized and accepted all terms and contents of the product's instructions, disclaimers and safety instructions. Users undertake to be responsible for their actions and all consequences. User undertakes to use the product only for legitimate purposes and agrees to these terms and any relevant policies or guidelines that the company may establish.
- ◆ In the process of using this product, please strictly abide by and execute the requirements including but not limited to the instructions and safety instructions. All personal injuries, accidents, property losses, legal disputes and other adverse events that cause conflicts of interest caused by violation of the safety instructions or irresistible factors shall be borne by users themselves, and the company shall not assume any responsibility.

please do not use wet hands to plug and unplug the power supply of the equipment.
 please be sure to use a regular brand power socket, and make sure the grounding is well grounded to prevent electric shock.
 please be sure to make regular safety checks on sockets and plugs to avoid potential electrical safety hazards caused by aging and short circuit.
 please be sure not to use the product in a humid or hot environment to ensure the safety of the product.
 please be sure to loosen bundle cords of the power cable before using, to avoid electromagnetic induction and heating, thus increasing the heat dissipation speed.
 please be sure not to use the product equipment or socket or other places in high or easy to fall, so as to avoid damage.
 before opening the socket power supply, please be sure to turn off the load power switch of the product equipment.
 the company reserves the rights to improve product, upgrade technology and change parameters without prior notice.