



←INSIZE→




**7302-8304**  
**Data output cable**  
**Manual**



产品介绍

**7302-8304:** Data output cable, suitable for **8304** series electronic balances

1. The data output cable RS232 port is connected to the electronic balance, and the USB end is connected to the computer for use, without the need to install any software or drivers
2. Open any Excel/WORD/TXT file, select a location with the cursor, and then press the transmission button of the data output cable or the "print" button of the electronic balance to transfer the weighing data to this location

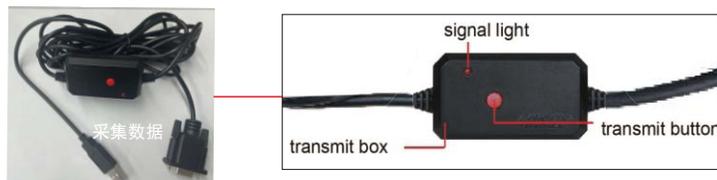

**7302-8304**  
Data output cable



**8304**  
Electronic balance

Explanation of electronic balance mode status transmission:

**8304-120/C:**

1. In C5-1 mode, data can only be transmitted by pressing the transmission button on the transmission box or the external foot switch (model: 7304-2, optional) on the transmission box, and cannot be transmitted by pressing the "print" button on the electronic balance
2. In C5-3 mode, only the "Print" button of the electronic balance can be pressed for transmission

**8304-220/C:**

1. In C5-1 or C5-3 mode, data can be transmitted by pressing the transmission button on the transmission box or by pressing the external foot switch (model: 7304-2, optional) on the transmission box, or by pressing the "print" button on the electronic balance

Note: The mode switching of the electronic balance can refer to the instruction manual of the electronic balance